


# SKOPED MICRO MICROSCOPE EYEPIECE SIZE FINDER




*Measure existing eyepiece. Printer needed.*

1. Print this page, be sure printer is set to print at 100%.
2. Verify the page printed correctly by first measuring the circle(s) diameter with a ruler, to ensure accuracy.
3. Measure the male end of the microscope eyepiece after you've removed it from your microscope, by putting it over the circle making sure the outside of the eyepiece lines up with the outside of the circle.
4. The circle that lines up closest, is the size of eyepiece you need to order. If it is 25mm or 28mm, these are special order pieces, so please reach out to our customer service email: [support@skopedmicro.com](mailto:support@skopedmicro.com) and we'll get you set up.

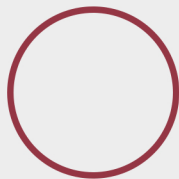


**23mm**

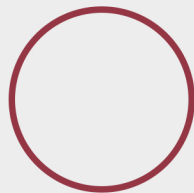


**25mm**

*\*Special order size*

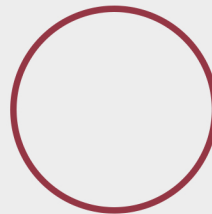

**28mm**

*\*Special order size*

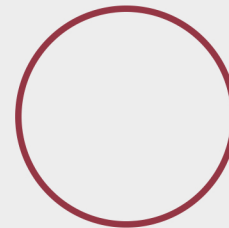

**30mm**



## Helpful Tips and Information

- There is a set screw that you may need to loosen to remove one of your eyepieces from your microscope.
- If you have multiple sizes of microscopes, we send a 30mm shim out with every 23mm eyepiece, to make it more universal. See our website for more information.
- Visit our YouTube Page @skopedmicro for more tutorials on how to find the correct size as well as how to use and get the most out of your Skoped Micro Setup.
- For any other questions call us at: **800-790-4480** or email us at: [support@skopedmicro.com](mailto:support@skopedmicro.com)

